


1. Configure the variable (plotting time) and click the button to specify the save file.

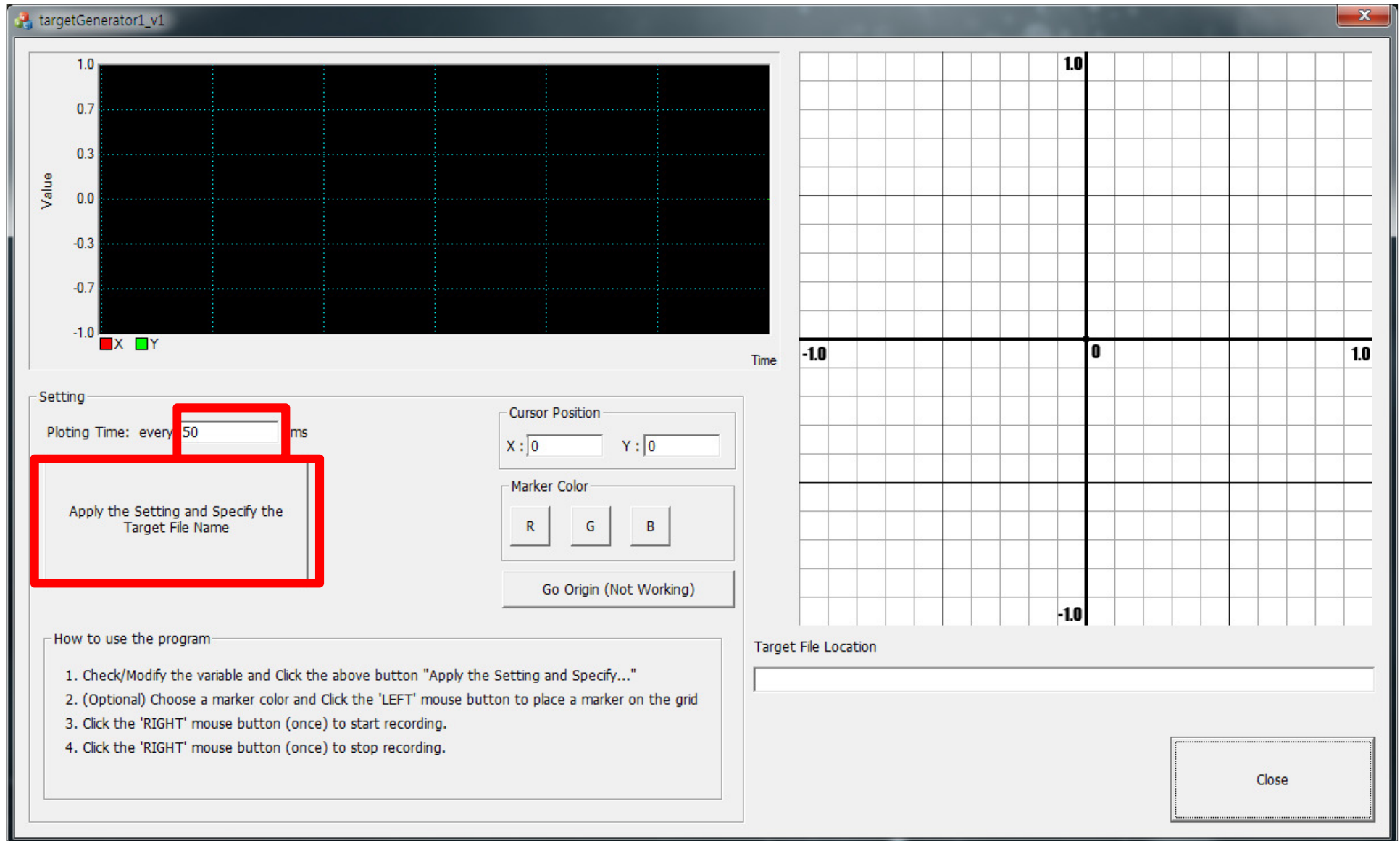

2. (Optional) If you want to leave a marker on the grid, choose the marker color (Red, Green, Blue) and LEFT Click on the grid.

The screenshot displays the 'targetGenerator1_v1' application window. It features a main grid with a coordinate system ranging from -1.0 to 1.0 on both axes. The grid is divided into four quadrants by a vertical line at x=0 and a horizontal line at y=0. The origin (0,0) is labeled. The axes are labeled 'Value' (vertical) and 'Time' (horizontal). The grid has a black background with a cyan dotted grid. A legend at the bottom left of the grid shows a red square for 'X' and a green square for 'Y'. Below the grid, there are several control panels:

- Setting:** A panel with a 'Ploting Time' field set to '50 ms' and a button labeled 'Apply the Setting and Specify the Target File Name'.
- Cursor Position:** A panel with 'X:' and 'Y:' input fields, both set to '0'.
- Marker Color:** A panel with three buttons labeled 'R', 'G', and 'B'. The 'B' button is highlighted with a red border.
- Go Origin (Not Working):** A button located below the Marker Color panel.
- How to use the program:** A text box containing four numbered instructions:
 1. Check/Modify the variable and Click the above button "Apply the Setting and Specify..."
 2. (Optional) Choose a marker color and Click the 'LEFT' mouse button to place a marker on the grid
 3. Click the 'RIGHT' mouse button (once) to start recording.
 4. Click the 'RIGHT' mouse button (once) to stop recording.
- Target File Location:** A text input field.
- Close:** A button in the bottom right corner.

The main grid shows several colored markers: a red marker at approximately (0.8, 0.3), a green marker at approximately (-0.5, -0.5), a blue marker at approximately (-0.8, -0.8), and another red marker at approximately (0.8, -0.8). The axes are labeled with '1.0' and '-1.0' at their respective ends.

3. If you click RIGHT mouse button **once**, then the recording starts. If you want to stop recording, click RIGHT mouse button **once again**.

